

Cat. # **WK551S**

For Research Use

TAKARA

GT-T551 Culture medium, 1L Bottle

Product Manual

v201602

Table of Contents

| | | |
|------|------------------------|---|
| I. | Description..... | 3 |
| II. | Component | 3 |
| III. | Storage | 3 |
| IV. | Note | 3 |
| V. | Application..... | 4 |
| VI. | References..... | 7 |
| VII. | Related Products | 7 |

I. Description

GT-T551 Culture medium is used for extended culture of human T lymphocytes as well as during T cell transduction by retroviral or lentiviral vectors. The medium contains only human serum albumin (pharmaceutical grade) and recombinant human insulin; no additional proteins or growth factors are added. GT-T551 Culture medium is tested to ensure it is sterile, free of endotoxins and mycoplasma, and promotes adequate lymphocyte proliferation activity. Correct pH and osmolality are also verified.

For T cell expansion, peripheral blood mononuclear cells (PBMCs) are typically used as starting material. Further T cell purification from PBMCs is not needed for T cell expansion. PBMCs can be separated by density gradient centrifugation from heparinized blood or by apheresis. Alternatively, it can be purchased from vendors.

II. Component

GT-T551 Culture medium, 1L Bottle 1,000 mL

III. Storage 2°C to 8°C**IV. Note**

GT-T551 Culture medium is supplemented with human serum albumin, human insulin, L-glutamine, and streptomycin, and it can be directly used for T cell culture. Although GT-T551 Culture medium is capable of supporting T cell proliferation without adding serum and plasma, cell growth may be enhanced by addition of these components in the medium. T cells can be cultured in GT-T551 Culture medium with cytokines (IL-2, IL-15, etc.) in an anti-CD3 antibody-stimulated state.¹⁻³⁾ During anti-CD3 antibody stimulation, the presence of RetroNectin® enhances the proliferation of T cells.⁴⁻⁸⁾ The expanded cell population shows a higher proportion of naïve T cells than anti-CD3 antibody stimulation alone (see page 6).

V. Application


GT-T551 Culture medium was used during T cell expansion with RetroNectin (Cat. #T100A/B) stimulation in a gas-permeable culture bag.

Note: The method of T cell expansion by RetroNectin stimulation requires a license from Takara Bio Inc. for uses other than research.

【 Method 】

- 1) Coat CultiLife™ 215 Culture bag with anti-CD3 monoclonal antibody and RetroNectin (Set-up; Day 0)
 - Prepare 30 mL of PBS containing anti-CD3 mAb (5 μ g/mL) and RetroNectin (25 μ g/mL).
 - Pour the solution into a CultiLife 215 Culture bag.
 - Incubate for 2 - 5 hours at 37°C in a 5% CO₂ incubator.
 - Remove the solution and gently rinse the inside of the bag with 30 mL of PBS.
 - Repeat the rinse step twice.

Note: On the last rinse, remove the wash solution just before adding PBMCs.
- 2) Add PBMCs
 - Suspend about 3×10^7 PBMCs in 30 - 50 mL of GT-T551 Culture medium containing human serum or plasma in a final concentration of 0.5 - 1.0%.
 - Pour the PBMC suspension into the anti-CD3 mAb-coated and RetroNectin-coated CultiLife 215 Culture bag prepared during Step 1).
 - Add GT-T551 Culture medium up to 200 - 300 mL and IL-2 (final concentration of 200 - 1,000 U/mL).
 - Culture at 37°C in a 5% CO₂ incubator.

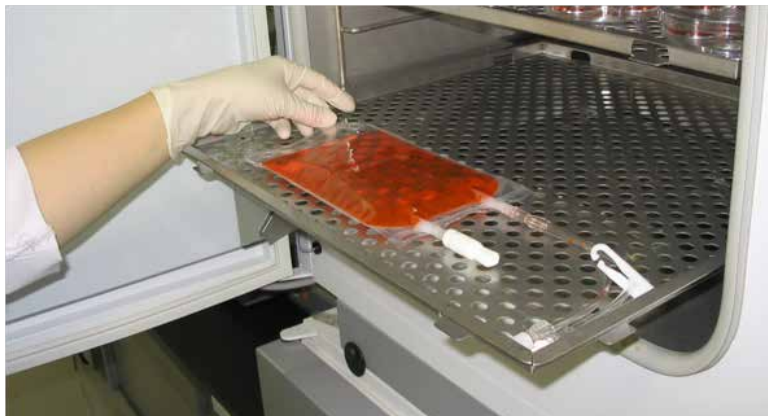


3) Passage at Day 4

- Collect all of the cell suspension from the CultiLife 215 Culture bag using the method described below and transfer to a new GT-T610 (CultiLife Eva) Culture bag (Cat. #FU0010).*

Note: Cells cultured in the RetroNectin-coated bag adhere to the bag.

Cells can be suspended by rocking the culture bag back and forth with both hands about 50 times as shown below.

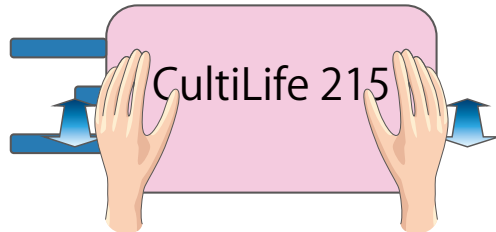

- Confirm by microscopy that no cells remain in the CultiLife 215 Culture bag.
- Add an appropriate volume of GT-T551 Culture medium and IL-2 (final concentration of 200 - 1,000 U/mL).

Note: Max volume of GT-T610 (CultiLife Eva) Culture bag is 1,000 mL.

- Culture at 37°C in a 5% CO₂ incubator.

* At this time, scale-up is possible by passaging the cells into more culture bags according to the rate of cell growth and the desired endpoint cell number.

4) Harvest cells

- Suspend cells by rocking the Culture bag about 20 - 30 times using the same method described in Step 3) for the CultiLife 215 Culture bag.
- Collect the cell suspension from the GT-T610 (CultiLife Eva) Culture bag.
- Wash cells by centrifugation, or if the culture volume is large, by using a cell washing apparatus such as the Cell Server.

* PBS or saline containing 0.1% human serum albumin can be used for the wash solution.

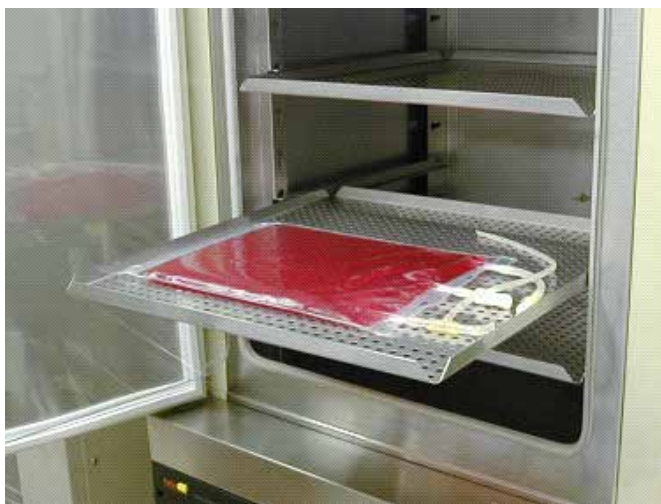


Table 1. Example of culture

| | Day 0 | Day 4 | Day 7 | Day 10 |
|----------------|-----------------------------------|---------------------------------------|----------------------------|---------|
| PBMC number | 3×10^7 cells | | | |
| Culture volume | 300 mL | 500 mL x 5 ^{*3} | 1,000 mL x 5 ^{*4} | |
| Culture area | 215 cm ² ^{*1} | 640 cm ² ^{*2} x 5 | | |
| | CultiLife 215 x 1 | GT-T610 x 5 | | Harvest |

*1: Culture area of one CultiLife 215 Culture bag is 215 cm².

*2: Culture area of one GT-T610 (CultiLife Eva) Culture bag is 640 cm².

*3: If the concentration of the cell suspension is less than 4×10^5 cells/mL at Day 4, it is recommended to passage on Day 5.

If passaging on Day 4, use a cell suspension concentration of at least 5×10^4 cells/mL.


*4: At this stage, passage using a cell suspension concentration of at least 5×10^5 cells/mL.

【 Example of T cell expansion using healthy donor PBMCs 】

Method

PBMCs and autologous plasma were prepared using 50 mL of blood from two healthy volunteers on the basis of informed consent. The effect of RetroNectin (RN) co-stimulation on T cell expansion was tested using the culture system above. Heat-inactivated plasma at a final concentration of 0.6% was added to the GT-T551 Culture medium at Day 0 and Day 4.

Results

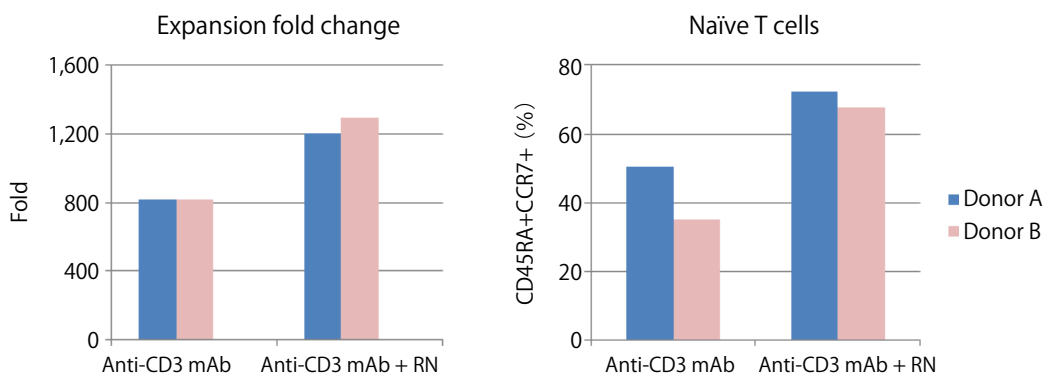

RetroNectin co-stimulation resulted in a dramatic increase in T cell expansion and the population contained a high proportion of naïve T cells.

VI. References

- 1) Wang, Y., *et al.* (2013) CIK cells from recurrent or refractory AML patients can be efficiently expanded in vitro and used for reduction of leukemic blasts in vivo. *Exp Hematol.* **41**(3): 241-252.
- 2) Ai, Y. Q., *et al.* (2014) The clinical effects of dendritic cell vaccines combined with cytokine-induced killer cells intraperitoneal injected on patients with malignant ascites. *Int J Clin Exp Med.* **7**(11): 4272-4281.
- 3) Dodo, K. *et al.* (2014) An efficient large-scale retroviral transduction method involving preloading the vector into a RetroNectin-coated bag with low-temperature shaking. *PLoS ONE.* **9**(1): e86275.
- 4) Yu, S. S., *et al.* (2008) In vivo persistence of genetically modified T cells generated ex vivo using the fibronectin CH296 stimulation method. *Cancer Gene Ther.* **15**(8): 508-516.
- 5) Ishikawa, T., *et al.* (2014) Phase I clinical trial of fibronectin CH296-stimulated T cell therapy in patients with advanced cancer. *PLoS ONE.* **9**(1): e83786.
- 6) Hosoi, H., *et al.* (2014) Stimulation through very late antigen-4 and -5 improves the multifunctionality and memory formation of CD8⁺ T cells. *Eur J Immunol.* **44**(6): 1747-1758.
- 7) Sakamoto, N., *et al.* (2015) Phase I clinical trial of autologous NK cell therapy using novel expansion method in patients with advanced digestive cancer. *J Transl Med.* **13**: 277.
- 8) Li, W., *et al.* (2015) Efficacy of RetroNectin-activated cytokine-induced killer cell therapy in metastatic brain tumor patients. *Oncol Res Treat.* **38**(4): 160-165.

VII. Related Products

RetroNectin® Recombinant Human Fibronectin Fragment (Cat. #T100A/B)
Anti-CD3 mAb GMP grade Anti-CD3 monoclonal antibody (Clone OKT3) (Cat. #T210)
CultiLife™ 215 Culture bag (Cat.#FU0005)
GT-T610 (CultiLife™ Eva) Culture bag (Cat. #FU0010)

RetroNectin is a registered trademark of TAKARA BIO INC.
CultiLife is a trademark of TAKARA BIO INC.

NOTE: This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc.

Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from TAKARA BIO INC.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.
